


96.96 Dynamic Array™ IFC

– Gene Expression

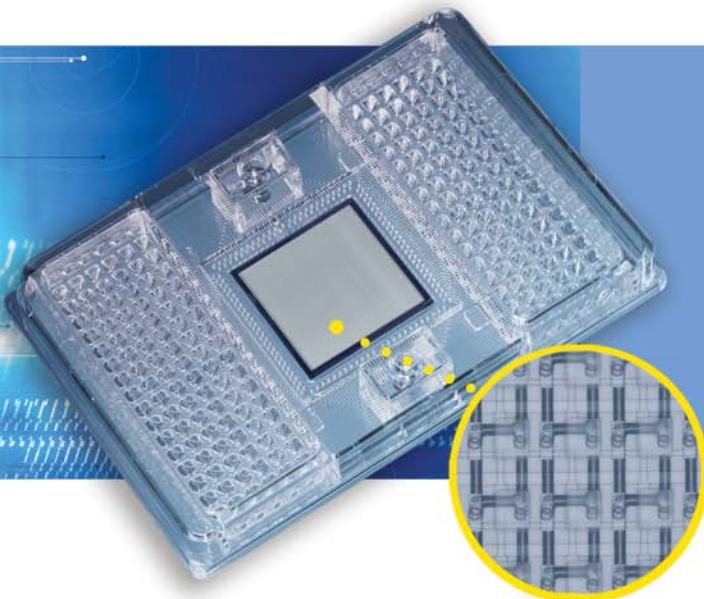

HIGH-THROUGHPUT MULTIPLEX PCR

The Fluidigm 96.96 Dynamic Array IFC – Gene Expression provides the flexibility of a microwell plate and the density of a microarray in one easy-to-use, consumable integrated fluidic circuit (IFC).


Key Benefits –

- Easy multiplexing of 96 primer-probe sets against 96 samples
- 9,216 individual data points per Dynamic Array IFC
- 192 liquid-transfer steps per 9,216 reactions, with complete setup flexibility

The New Standard in High Throughput Profiling

Fluidigm Dynamic Array IFCs radically reduce the cost per data point and time to results while radically raising the bar for parallel throughput. The chart below shows parameters to complete a study of 2,000 samples against 96 genes using 384-well plates compared to 96.96 Dynamic Array IFCs:

	384-WELL MICROPLATES	96.96 DYNAMIC ARRAY IFCs
TOTAL RUNS	500	21
REACTIONS PER RUN	384	9,216
TOTAL LIQUID-TRANSFER STEPS	384,000	4,032
TOTAL MASTER MIX	960 ml	5.1 ml

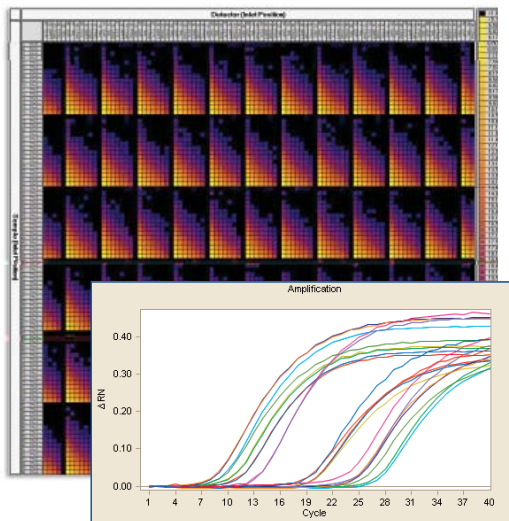

Gene expression results may be viewed as real-time curves and as a heat map showing 9,216 reactions per run.

The Power of Microfluidics

With a Dynamic Array IFC, high-throughput multiplexing is easy because the microfluidic architecture does the work of combining samples and primer-probe sets into 9,216 PCR reactions. That's twenty four-fold more data than is produced by a 384-well plate. This radical advance in experiment density is fully leveraged through a hardware/software system that automates setup and data analysis.

96.96 Dynamic Array IFC – Gene Expression

Specifications

PARAMETER


Quantitative resolution	2-fold difference in starting copy with 99.7 % confidence and 6-log of dynamic range
Chip format	SBS Compatible (128 mm x 85 mm x 14 mm)
Inlet spacing on input frame	4.5 mm pitch
Liquid transfer steps	192
Primer-probe inlets	96
Sample inlets	96
Reaction chambers	9,216
Reaction volume	6.7 nl
Instrument compatibility	BioMark Real-Time PCR System, IFC Controller HX

Fluidigm System for Genetic Analysis

- **Dynamic Array™ IFCs**
Consumable IFCs for high-throughput gene expression analysis and SNP genotyping.
- **Digital Array™ IFCs**
Consumable IFCs for digital PCR.
- **IFC Controller**
Integrated hardware/software for loading IFCs.
- **FC1™ Cyclizer**
Integrated hardware/software for thermal cycling of IFCs.
- **EP1™ Reader | Real-Time PCR System**
Integrated hardware/software for detection of fluorescent signal within IFCs.
- **Software Suite**
Analysis software for gene expression analysis, SNP genotyping, and digital PCR.
- **Service Plans**
Hardware service and software maintenance plans.

Work Flow

- 1 Prime**
Prime the Dynamic Array IFC to close the interface valves, preventing premature mixing of samples and assays.
- 2 Transfer**
Pipette samples, premixed with master mix, into separate sample inlets and the primer-probe sets into separate primer-probe inlets on the IFC.
- 3 Load**
Place the Dynamic Array IFC on the IFC controller, and use the software interface to pressure load the assay components into reaction chambers. Assay components are automatically combined on-chip.
- 4 Run**
Place the Dynamic Array IFC on the BioMark Real-Time PCR System for thermal cycling and fluorescence detection.
- 5 Analyze**
Use real-time qPCR Analysis software to view and to interact with amplification curves, color-coded heat maps, and C_t data for the run.

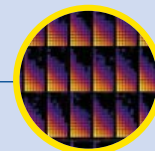

For Use with Gold-Standard PCR Assays

The BioMark system runs licensed 5' nuclease assays, so it integrates easily into established workflow. The footprint of the Dynamic Array IFC and spacing of fluid inlets comply with SBS* standards, so the laboratory may continue to use existing liquid-handling equipment. Fluidigm has adopted SBS standards for all of its systems, ensuring compatibility of BioMark instrumentation with higher density IFCs in future releases.

Fluidigm®

Corporate Headquarters

Fluidigm Corporation
7000 Shoreline Court, Suite 100
South San Francisco, CA 94080 USA
Toll-free: 1.866.FLUIDLINE
Fax: 650.871.7152
www.fluidigm.com

Sales

North America
650.266.6170 | biomark@fluidigm.com
Europe/EMEA
+33 1 60 92 42 40 | biomark@fluidigm.com
Japan/Korea
+81 3 3555 2351 | biomarkasia@fluidigm.com
Asia
biomarkasia@fluidigm.com

* The Society for Biomolecular Sciences

© Fluidigm Corporation. All rights reserved.
Fluidigm, the Fluidigm logo, BioMark, EP1, FC1, Dynamic Array, Digital Array, and FLUIDLINE are trademarks or registered trademarks of Fluidigm Corporation in the U.S. and/or other countries. All other trademarks are the property of their respective owners.
Fluidigm recommends that you only purchase TaqMan® dual-labeled probes and/or other licensed PCR assay reagents from authorized sources.